

Food Packaging Residual Solvents

Column: **MEGA-1701 - 0.15mm, 1.00µm, 10m**

Catalog Code: S-1701-015-100-10


Conditions:

Injection: Split 250°C, 0.2µL with micro-syringe,
200mL/min Split Flow.

Detector: FID 250°C.

Oven Temp.: 35°C (0.5 min), 13°C/min, 110°C.

Carrier Gas: Helium, 140 kPa.

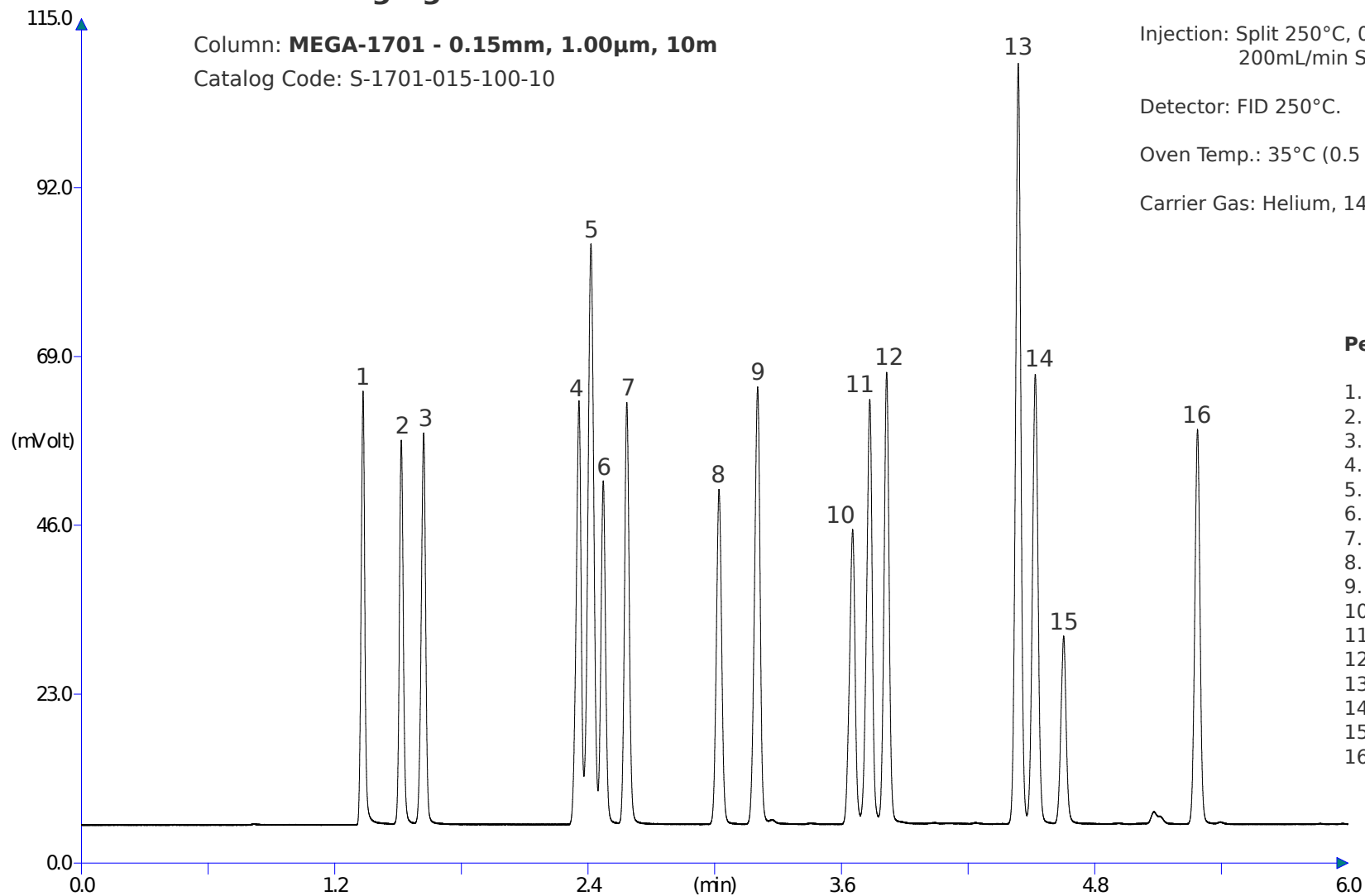

Peaks:

1. Ethanol
2. Acetone
3. 2-propanol
4. 1-propanol
5. Cyclohexane
6. Ethyl acetate
7. Methyl ethyl ketone
8. Isopropyl acetate
9. Isobutanol
10. Methoxy propanol
11. 1-butanol
12. Propyl acetate
13. Toluene
14. Isobutyl methyl ketone
15. Ethoxy propanol
16. Butyl acetate

